
Leica EM ACE200 Sputter Coating: Quick Start Guide for Technical Staff

V1: S. Herrmann, 12/20/2013

~~V371T-V374~~ | ~~July 2017~~ March 2018

Introduction

The **Leica EM ACE200** coating system is used for precise coating of samples for subsequent examination with a scanning electron microscope (SEM). Samples are metal coated using the sputtering method where argon plasma erodes a target material. Carbon coating is achieved by carbon thread evaporation. Any material can be processed as long as it is not sensitive to vacuum, argon plasma, or the heat generated during carbon coating. Sputter coating prevents charging the specimen with an electron beam in the SEM high-voltage mode. Metal coatings are also useful for increasing signal-to-noise ratio. The ACE200 can be configured as a sputter coater, carbon coater, or both, where the two processes can be switched over from one to the other.

With the Hitachi TM3000 SEM, sputter coating is not required for all samples. Images up to a magnification of ~~×4000x~~ are of high quality without any sputter coating in **Charge-up Reduction Mode (Settings > Observation Mode)**.

For detailed observations in magnifications greater than ~~×4000x~~ and for very delicate fossils, sputter coating before using the SEM can be used to potentially receive higher quality images.

Apparatus and Materials

The Leica EM ACE200 coating system includes the following main functional units (**Figure 1**):

- Vacuum chamber
- Touch screen control panel
- Sample stage with 18 positions for 1/2” SEM stubs
- Removable internal shielding, shutter, and door
- Carbon source and sputter targets
- Vacuum pump
- Sample stage



Figure 1. Leica EM ACE200 Sputter Coater

Targets

The following targets are used for sputter coating:

- Gold
- Gold-Palladium
- The instrument can use other metal targets but these are not in use at this time

Argon gas supply

The working gas (argon) must be supplied under a pressure of ~ 500 mbar ± 100 mbar (5.6—8.4 psi above atmosphere). The gas should be at least 99.99% pure.

Caution! Do not use a standard two-stage regulator for the sputter coater unless it is equipped with a low-pressure delivery-side gauge. Most regulators don't effectively pick up pressure until 10 or even 15 psi or higher. It may be necessary to put a second low-pressure-range regulator downline from the main line regulator. Make sure you have the right tool for the job!

Sputter coating

Sputter coating is performed using ionized argon to create a plasma. The argon ions are accelerated by high voltage and directed toward the source via a magnet where they collide with the target and displace surface atoms. Due to this collision the surface atoms are directed toward the area below the target and coat the sample. This coating process can be set to directional or diffuse (which provides more even coating on a bigger surface and is better for fissured samples), depending on the process pressure. This also influences the coating rate (diffuse means slower rate) and the grain size (directional means finer grains).

Loading the sputter target

There is an ACE200 youtube video that shows the loading of the target at <http://www.youtube.com/watch?v=Kd1lyKIDT8>.

The procedure is as follows:

1. Wear gloves and make sure the instrument is fully vented.
2. Open the source cover and unplug the connectors.
3. Unscrew the 2 sputter head screws to remove the flange.
4. Remove the bayonet ring by turning and insert the sputter target as shown in Figure 2, below.
5. Fix the target by gently tightening the bayonet ring; **tighten by hand only**.
6. Replace the sputter head and gently tighten the fastening screws.
7. Connect the cables and close the cover.

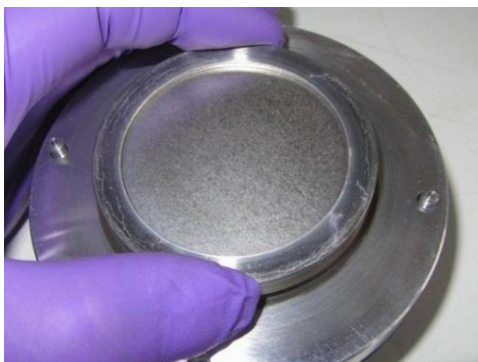


Figure 2. Inserting the sputter target.

Setting up a Sputtering Process

Sputter Parameters

Target material: select from a drop-down list; the instrument automatically chooses parameter settings.

Directional mode: intended for relatively flat samples; base vacuum is 4×10^{-2} mbar for this mode.

Diffuse mode: intended for more topographic samples; base vacuum is 8×10^{-2} mbar for this mode.

Sputter current/purge cycles: automatically set when target is selected. Change under the **Setup** button.

Process termination time: select the **Clock** icon then use the +/- buttons to select a time in seconds.

Rotation speed: press the **Rotation** button, check **Rotation On**, and use the +/- buttons to set the speed.

The sputtering procedure is summarized as follows:

1. Turn on the main power supply to the instrument and make sure the chamber door is closed, the cover is closed, and the sputter head is connected (see **Loading the Sputter Target**, above).
2. Select the **Sputtering** button.
3. Press the **Pump** button to start pump-down.
4. During pump-down, set the process parameters:
 - a. Target material (last material used is presented by default)
 - b. Mode (**Directional** or **Diffuse**) button
 - c. Sputter current/purge cycles (preset when target was selected)
 - d. Process termination time (e.g., 12 seconds)
 - e. Rotation speed (e.g., 1)
5. Press the **Start** button (which then becomes the **Stop** button).

Note: if the vacuum cannot stabilize, make sure the argon supply valve is open and pressure is on the line.

Sputter Process Start

The coating process can be started at any time (system evacuated or not) as long as the door and source cover are closed and the source is connected. The **Start** button will turn into a **Stop** button when the system starts the coating cycle.

Once the **Start** button is activated the coater automatically runs the complete coating cycle. The system either stays under vacuum or vents automatically (if **Vent after process** is activated).

Pressing the **Stop** button terminates the process after confirmation, regardless of the step of the process. The system either stays under vacuum or vents automatically.

After setting coating parameters and pressing **Start** the system performs the following steps automatically:

1. Pump until base vacuum is reached.
2. Stabilize the plasma.
3. Pre-sputter, if target requires it (to clean oxidation from the target and to enable a stable sputter rate).
4. Start the sputter processing by opening the shutter and starting rotation (if activated).
5. Terminate sputtering by time.
6. Close the shutter.
7. Display results of the process.
8. Vent or stay under vacuum, per settings.

Carbon thread coating

The carbon coating process is carried out by evaporating a carbon thread. It is possible to coat using short pulses of 150 milliseconds or evaporating the thread completely with maximum power, a so-called flash.

Loading a carbon thread

The carbon thread can be loaded as a single thread or as a double thread. Thin layers from 1 to at least 20 nm can be achieved (there is variance in threads).

To minimize carbon thread waste, when loading a double thread, cut a piece of thread twice as long as the width of the black door frames of the coater. Fold the thread into half and load it as follows:

0. **Before beginning**, prepare the following on a clean surface: gloves, carbon head, Torx TX 10 key, brush, carbon thread.
1. Make sure the instrument is vented. Wear gloves.
2. Open the source cover and unplug connectors.
3. Unscrew the 2 evaporation head screws to remove the flange.
4. Loosen all 5 clamp screws with the Torx key.
5. Remove any carbon residue using the brush.
6. Loop the carbon thread around the first clamp; pull both ends gently to the left while tightening the screw.
7. Wind the thread around the other clamps, taking care that the thread slides into the clamping groove.
8. Pull the thread gently and tighten the last clamp.
9. Tighten the 3 remaining screws and trim thread on both ends as shown in Figure 3, below.
10. Replace the head, gently tighten the fastening screws, connect the cables, and close the cover.

Also note that there is an ACE200 youtube video where the loading of the thread can be seen at <http://www.youtube.com/watch?v=yKd1lyKIDT8>.

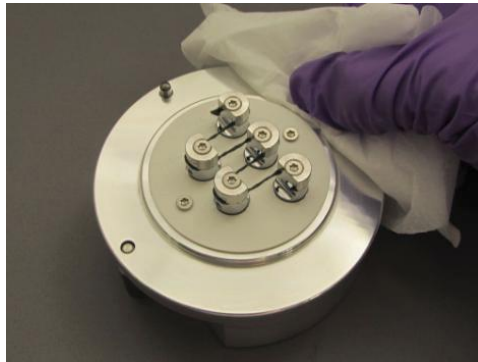


Figure 3. Threading the Carbon Thread

Processing Carbon Thread Coating

After setting coating parameters and pressing **Start** the system performs the following steps:

1. Pump until base vacuum is reached.
2. During pump-down, the following parameters will be set:
 - a. Target thread will be set to 2 if double-thread (automatic)
 - b. Process pulse/flash: use the +/- buttons to set number of pulses or flashes
3. Check availability of at least one carbon thread section.
4. Outgas (preheat) the first carbon thread for 45 seconds.
5. Open shutter and start rotation.
6. Pulse/flash according to the settings.
7. Close the shutter.
8. Display results of the process (number of pulses/flashes, availability of further threads).
9. Vent or stay under vacuum.

Switching between Processes

To switch between carbon thread and sputter coating, exchange the source (sputter or carbon thread) and choose the process on the main screen of the instrument. The instrument automatically detects if the correct source is installed and gives an error message if it does not match the selected process.

Safety

All electronic components are protected by covers (door, source cover, housing). The door and source covers are equipped with sensors that cut off power when they are opened. Software interlocks cut off power when a malfunction is detected (e.g., vacuum leak, short circuit in the source head, missing contact in the system).

Overheat protection: at 65°C the process is discontinued until the temperature drops to 45°C, then the process continues.

In case of a sudden vacuum breakdown, the instrument switches off automatically to protect pump and electric parts.

If the coating system is damaged or malfunctions, all use of the system should be suspended until the malfunction or damage has been corrected.

Do not operate the coating system unless all covers are properly in place – there is danger of electric shock and burns from high-temperature components.

Maintenance (Technical Staff)

For all maintenance the system has to be switched off.

Wear gloves when working on components in the vacuum area and for adjustment work.

Change the vacuum pump oil once per year.

Tighten hinges on the cover using a 3 mm Allen key if the cover closes too quickly or is loose.

Clean or replace shutter when a peeling or thick coating is visible, if contamination is observed on SEM, or vacuum pump-down time increases significantly (vacuum should reach 5 mbar after 30 sec pump-down).